

SLT4/SLT5 Series Adjustable Height Science Lab Table Instructions

The SLT4/SLT5 Series has been designed with the capability to be used with both our standard non-book compartment style table or with book compartments (purchased separately).

Note: *Extra apron marked BC is for use with book compartments only.*
For optional book compartment units, proceed to pg. 2 for book compartment installation instructions which should be done prior to leg installation.

Follow the below instructions for installation of adjustable legs.

Tools Required: Power screw driver with Philips head

Step 1: Place the tabletop upside-down on the floor with apron side up.

Note: Floor should be clear of any objects or dirt in order not to damage the tabletop.

Step 2: Carefully remove each of the four height adjustable legs from second carton and align holes on leg mounting plate with pre-drilled holes at the four corners of the underside of the table.

Step 3: Using a screw driver, carefully install each of the four legs with the provided screws (6 screws per leg) and tightly secure each of the screws.

Step 4: Attach a glide to the bottom of each leg by inserting the threaded end of the glide into the threaded hole on the bottom of the leg.

Step 5: Turn the table over with the tabletop facing up. Place the table in desired location. Adjust the glides to assure that table is stable.

Book compartment installation instructions.


Removing Standard Apron Panel

Place the tabletop with apron side up on a clean surface.

Note: Surface should be clear of any objects or dirt to avoid any damage to tabletop.

Using a screwdriver, begin by removing one front apron panel from tabletop in the following manner:

- Remove the screws that secure the apron panel to the tabletop (see figure A)
- Remove the screws that secure the front apron panel to the brackets. (see figure B)
- Remove the screws from the two apron panel reinforcement bars that attach to underside of table. (see figure C)

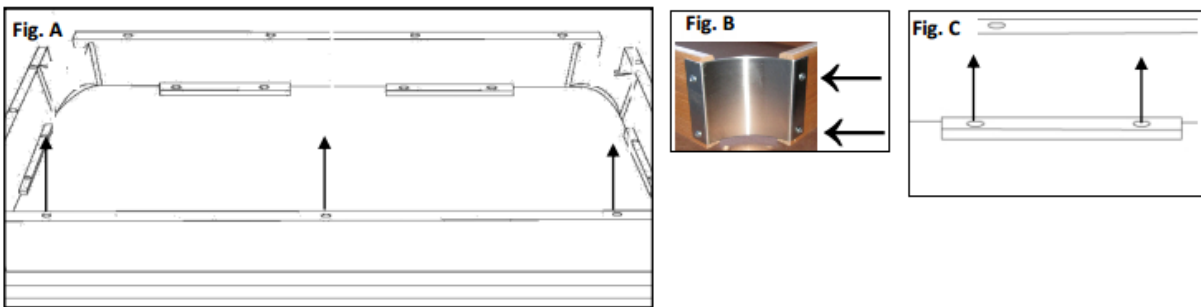

Installing Book Compartment Apron Panel


Remove apron panel marked **BC** from tabletop carton and install onto tabletop in the following manner:

- Place the BC apron panel in the location where the original apron was located and attach the brackets with two screws on each side.
- Further secure apron by inserting the screws from the front panel into the tabletop.

Installing Book Compartments

Place book compartments upside down on the tabletop and insert the bottom firmly into the grooves on the apron panel (see figure D)

Using a Philips screwdriver, fasten twelve screws (for each compartment) into the holes of the book compartment reinforcement bars. (See figure E)

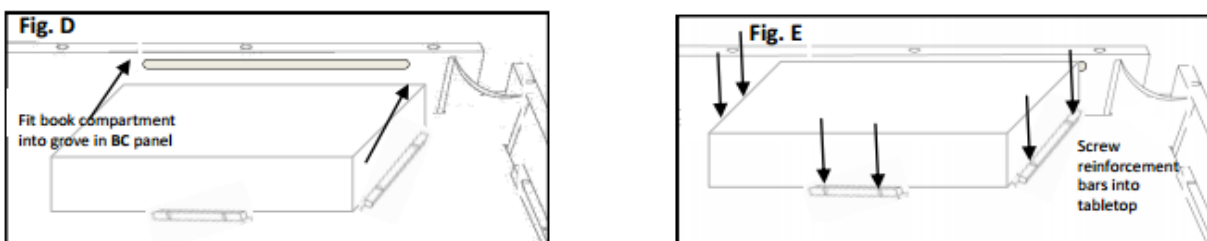

Once book compartment installation has been completed, proceed to install legs by following above instructions.

For optional caster installation, unscrew glides and tightly screw casters to bottom of legs.

CAUTION: Double check all screws to be tightened securely before use. Check all screws periodically.

Instructions for table height adjustment

1. For incline use, stand at one side of table, place feet on extended portion of glides and lift to desired height. Legs will automatically lock in place. See Ill. 1. For full table adjustment, repeat this at other side of table.
2. To lower table, stand at one side of table, place feet on extended portion of glides and lift to the highest height position then pull upwards to release. Gently push tabletop and lower to bottom-most position. Repeat this at other side of table.



Dear Customer,

Thank you for your purchase. Although extreme care goes into manufacturing & packaging of this product, mistakes are possible. If you encounter any missing parts or difficulty in assembly, please contact our customer service **Toll Free at 1-800-261-4112** or info@nationalpublicseating.com for courteous & immediate resolution to your problem.